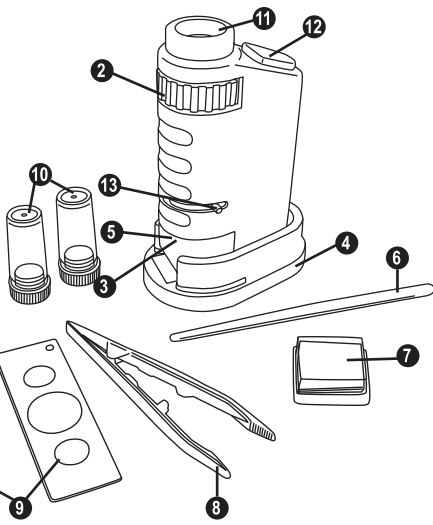
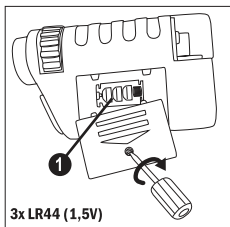
Mini-Mikroskop-Set

19444

(D)

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(GB)

With 'der echte Scout' mini microscope, you can view tiny objects close up.

Inserting the batteries

Open the battery compartment (1) on the side of the microscope and insert the batteries 3x LR44 (1,5V) (Batteries included.) as shown.

Then put the battery compartment cover back and screw it in place.

Important! Remove rechargeable batteries from the item before recharging! Recharge rechargeable batteries only under adult supervision! Non-rechargeable batteries are not to be recharged. Different types of batteries or new and used batteries are not to be mixed. Only batteries of the same or equivalent types as recommended are to be used. Batteries are to be inserted with the correct polarity (pay attention to the +/- pole). Exhausted batteries are to be removed from the item. The supply terminals are not to be short-circuited. Remove all the batteries when the item is not in use for a long period. Do not throw away old devices and used batteries in the household garbage. Disposal Instruction: Please only place discharged batteries into the collection boxes at the store or the municipal disposal site.

This toy includes test batteries. Please change against new batteries, if the batteries should be discharged and the function of the toy is affected.

Warning: Dispose of used batteries immediately. Keep new and used batteries away from children. If you think batteries might have been swallowed or placed inside any part of the body, seek immediate medical attention.

Function

A microscope is a device with which you can view tiny objects or capture images of these under

powerful magnification. The objects usually viewed and their structures are so small that they cannot normally be seen with the unaided human eye. Microscopes are important tools used by researchers in biology, medicine and the material sciences.

This mini microscope, with its stand (4) and accessories - including a stirring rod (6), a cleaning cloth (7), tweezers (8), two slides (9) and two storage containers (10) - allows you to see objects in 40x magnification.

Use

Place the microscope on the stand (4). Press the on/off button (12) on the microscope until the LED light (5) of the viewer lights up. Position what you want to view between the slides (9). Now place your eye on the eyepiece (11).

The lens separates light rays so that an optically enlarged intermediate image of the object on the slide (9) is produced on the viewer. The length of barrel of the microscope can be adjusted (13) using the focus adjustment (2) of the lens, so that, like a magnifying glass, the image of the object is enlarged. With the help of the tweezers (8), the stirring rod (6), the two storage containers (10), the two slides and the cleaning cloth (7) you can set up your own mini laboratory.

Care and warnings

Use the cleaning cloth to keep lenses and slides clean. Never use your fingers to clean the lenses or slides as this can result in scratching. Never force the microscope as although it and its accessories are robust, they can be damaged or broken.

Warning. Do not aim at eyes or face. Warning. Not suitable for children under 36 months. Small parts. Choking hazard. Do not directly look at the sun! Subject to technical change and change of color. Please keep this information for any possible correspondence. Warning: For safety reasons, please remove all packaging, wires, rubber bands, tags or other fastenings before handing this toy to a child.

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Happy People GmbH & Co. KG